

Use Instructions

1. Unwrap the enclosed meshes
2. Based on your food viscosity, decide which mesh hole size will best help remove food clumps
3. Do not bend the meshes! Simply drop your chosen mesh inside the strainer and remove by pushing up from the bottom
4. The strainer is designed to fit/attach on the back left or right side of your Droso-Filler
(See photos on our website: geneseesci.com)
5. Before using, heat the Droso-Strainer and mesh as you would the Droso-Filler. Both are necessary to help prevent food clumping
6. Set Droso-Filler upright
7. The strainer and meshes are made of durable stainless steel. They can be immersed and cleaned with soap and water using a soft non-scouring brush or nylon sponge. Both can be autoclaved



Customer Service

If you need assistance, please contact us at **858.536.8044** or email at support@geneseesci.com.